sunscreen

UNCHAINED
LABS

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
The screenshot displays the Sunny Suite software interface for an mRNA-LNP Demo-1 experiment. The interface is divided into several sections:

- Protocols:** A sidebar on the left lists various protocols, with 'mRNA-LNP Demo-1' selected.
- Configuration:** A central panel shows the experiment's description, apparatus (Input 1, Input 2, Collection), and parameters (Configuration: 1.000meters).
- Execution Time:** A large digital display shows '00:00:24'.
- Image:** A central image shows the Sunscreen hardware with a 'Configuration' overlay.
- Status:** A section below the image indicates 'Experiment 01/06: Preparation' and includes 'Start' and 'Stop' buttons.
- Execution Control:** A section below the status area includes radio buttons for 'Run all rows' and 'Run from row n'.
- Output:** A log on the right side shows a detailed timeline of the experiment's progress, including steps like 'Sunscreen Experimental Protocol Starting', 'Input 1 well plate definition assigned', 'Input 2 well plate definition assigned', 'Collection well plate definition assigned', 'Running all experiments', 'System preparation', 'Preparation', 'Dilution not used', 'System preparation', 'Filling system pump', 'Nucleoside filling', 'System pump: State set to Filling', 'System preparation', 'Filling system pump', 'System pump: Fill at rate 8000 µl/min', 'Nucleoside filling', 'System pump: State set to Filling', 'Input preparation of mRNA Acetate', 'Input preparation of SMA-102 Ethanol', 'Collection preparation', and 'System Pump: State Filling'.
- Manual Control:** A bottom section shows 'Current Readings' for 'Input 1', 'Input 2', and 'Collection', each with 'Pressure (System)' and 'Pressure (Sample)' values (e.g., 0.3 bar) and 'Bury'/'Stop' buttons.

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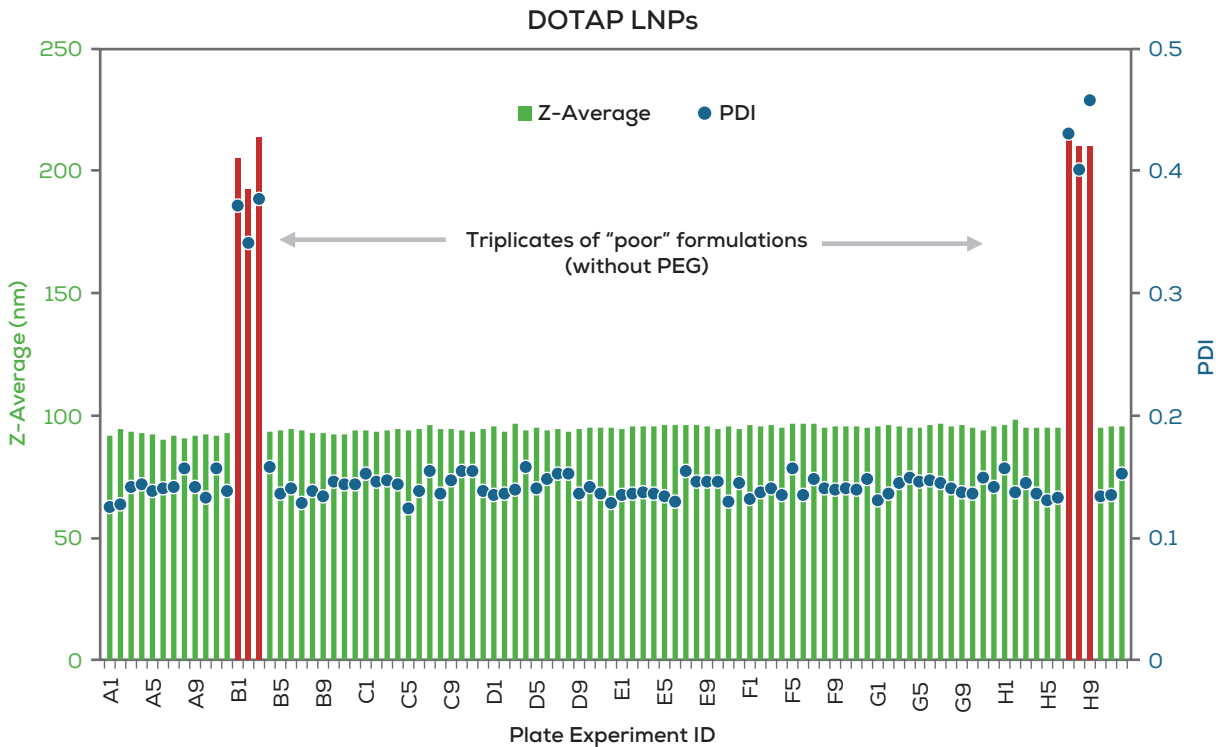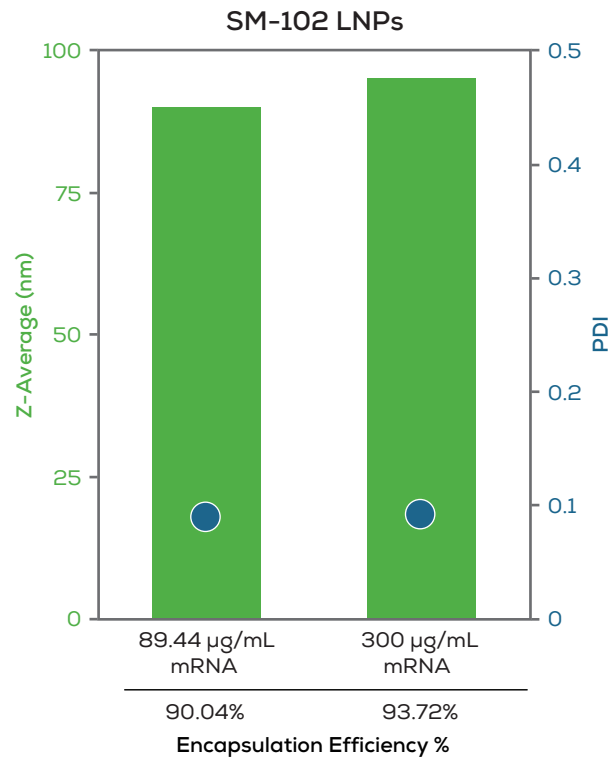
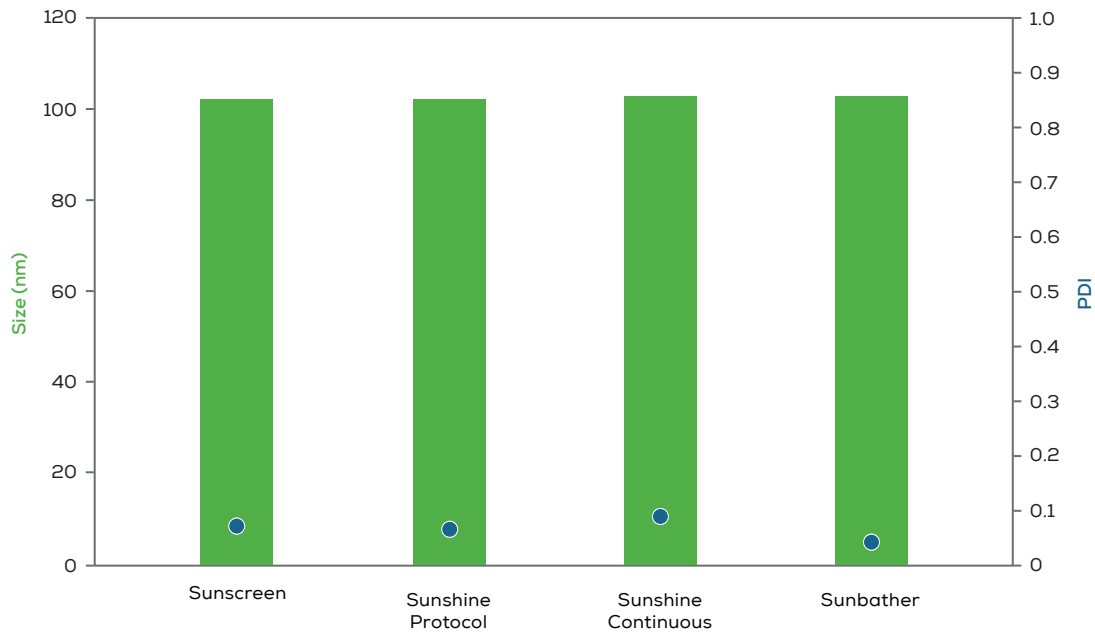
© 2024 Unchained Labs. All rights reserved. The Unchained Labs logo, Stunner, Sunscreen and the Sunscreen logo are trademarks and/or registered trademarks of Unchained Labs. All other brands or product names are trademarks owned by their respective organizations.