sunshine

UNCHAINED  
LABS

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The screenshot displays the Sunny Suite software interface for configuring and running a protocol. The main window is titled "Sunny Suite Version 1.0".

**Left Panel (Protocols):** Lists protocol families including "Sunshine Development", "Sunshine Prime", and "Sunshine System Initialize". The selected protocol is "BK Timing Test Sunshine - 22-SEP-23-1".

**Configuration Panel:** Shows the protocol description and apparatus. The apparatus table is as follows:

Function	Device
Automated sample valve	ASV
Input 1 pump	Aqueous
Input 2 pump	Ethanol
Dilution pump	Dilution
Automated collector	Automated Collector

**Execution Panel:** Shows the execution time as 00:04:14, an image of the system labeled "Configuration A", and the status "Sunshine protocol: System Preparation". The execution control section indicates "Experiment is running" with "Start" and "Stop" buttons. Below this, there are radio buttons for "Run all Experiments" and "Run from Experiment".

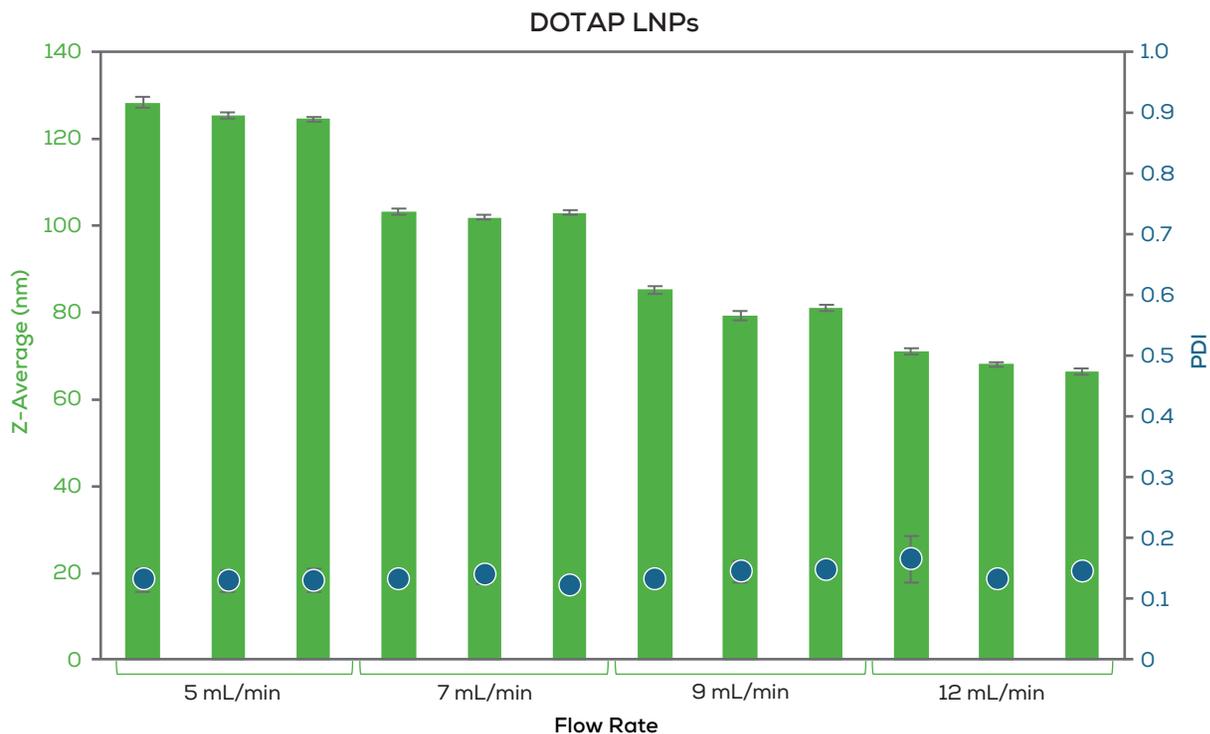
**Output Panel:** Displays a log of execution events, including "Sunshine protocol VQ 1.0.0 Test Automated system for mixing and fractionation of lipid nanoparticle formulations", "Sunshine protocol Starting", "Sunshine protocol Automated sample valve. Firmware version (1.14) is supported", "Sunshine protocol Automated collection Stop", "Sunshine protocol Lib All pumps fitted with Red Springs", "Sunshine protocol Lib Collector rack description selected 2 ml Rack, 35 Vials, 5x7 (7200130) matches imported rack description", "Sunshine protocol Running all experiments from selected experiment", "Sunshine protocol A System Preparation", "Sunshine protocol A-1 Collection system preparation", "Sunshine protocol A-1.1 Collection Vial to wash position", "Sunshine protocol Automated collection Positioning at Wash position", "Sunshine protocol A-1.2 Collection valve to COLLECT", and "Sunshine protocol Automated collector Setting valve to Collect".

**Bottom Panel (Manual Control):** Shows the status of various components: Automated Collector (Idle), ASV (Idle), Aqueous (Idle), Ethanol (Idle), Dilution (Idle), and V Quad Pump 4 (Not initialized). Each component has a "Current Readings" section with "Valve" and "Pressure" values. For example, the Aqueous pump shows a pressure of 0.3 bar.

**Bottom Right:** Displays system information: "RAM: 322MB" and "CPU: 4%".

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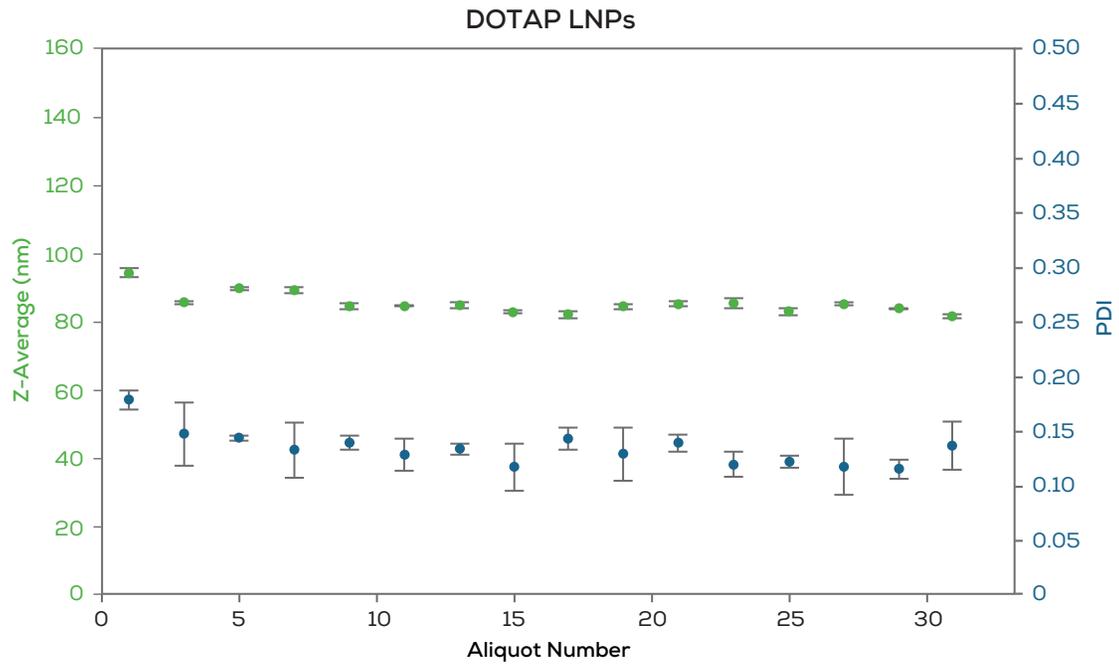
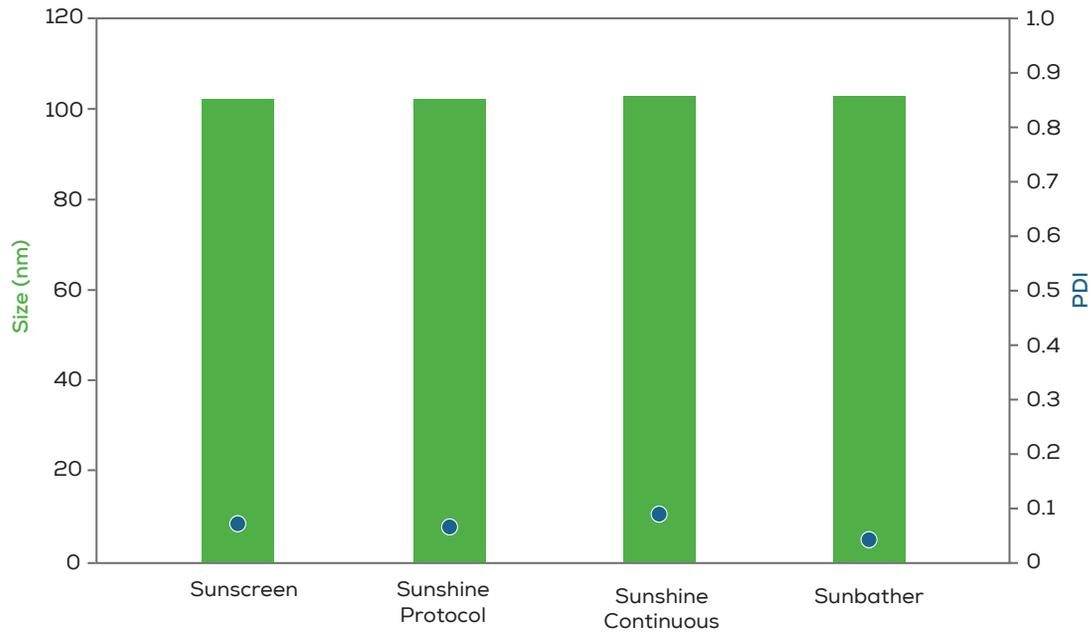
© 2024 Unchained Labs. All rights reserved. The Unchained Labs logo, Stunner, Sunscreen, Sunshine and the Sunshine logo are trademarks and/or registered trademarks of Unchained Labs. All other brands or product names are trademarks owned by their respective organizations.

Rev B